

Sanotex® Environmental Surface Wipes

Compatible with your preferred chemistry

These low-lint, disposable microdenier wipes pull from a reusable dispenser with a spring loaded lid to offer quick, reliable, cost-effective cleaning. Constructed of ultrafine meltblown fibers with a core of stronger spun-bonded fibers for optimal strength, Sanotex are compatible with peroxides, peroxide/peroxyacetic acid, phenolics, and quaternary ammonium-based formulations, and deliver chemicals to surfaces efficiently.



Features

Cost-effective, 100% synthetic, and disposable

Self-closing, spring-loaded lid keeps wipes wet longer

Measured chemical usage and flexible chemical choice

Compatible with bleach, quat, and peroxide-based disinfectants

Part No.	Description	Size	Packaging
85337	Sanotex Environmental Surface Wipes	10" x 12" (25 x 30.5 cm)	60/roll; 30 rolls/case
85338	Sanotex Environmental Surface Wipes	10" x 12" (25 x 30.5 cm)	40/roll; 8 rolls/case
85339	Sanotex Environmental Surface Wipes	6" x 8" (15.2 x 20.3 cm)	150/roll; 6 rolls/case
85340	Sanotex Environmental Surface Wipes	10" x 12" (25 x 30.5 cm)	60/roll; 6 rolls/case
85347	Sanotex Environmental Surface Wipes	10" x 12" (25 x 30.5 cm)	40/roll; 48 rolls/case
85357	Sanotex Environmental Surface Wipes	6" x 8" (152 x 203 cm)	150/roll; 30 rolls/case
85341	Sanotex self-closing dispensing canister	4.75" x 8" (120 x 200 cm)	6/case

How to use Sanotex® Wipes



< Insert a dry roll of Sanotex wipes into the canister



< Carefully pour chemical solution over center of wipe roll in canister. Allow five minutes before use to assure proper saturation of wipes.



< From the center of the roll, pull a wipe up and carefully thread through opening in canister lid (approximately one inch). Replace lid, screwing on tightly.



< To dispense, open the hinged lid and pull the wipe through the opening at an angle. The spring-loaded lid will reclose automatically.